FRANK-PTI

QUALITY TESTING INSTRUMENTS

DE/EN

40690 ProbeNet

Erstellt von: / Issued by: FRANK-PTI

Version: Version 1.0

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2. ENGLISH

2.1. General

This guide serves as a training material and makes it easier to operate the equipment properly and safely.

A professional operation and careful maintenance are essential for obtaining the reliability and operational capability of the machine. The safety of all persons that come into contact with the machine depends on the mastery of the machine.

Therefore:

Read this manual before using the machine for the first time!

The manual makes it possible to be acquainted with the machine and its possible uses prior to using it for the first time and, to use it for its intended purpose. It contains important information which ensures a functional, economic and safe operation of the machine.

The operating manual is not only an indispensable introduction for the operator that will be taught, but contains useful tips. It contains instructions and suggestions for experienced workers. It is a necessary reference book for all users of the machine. By reading the operating instructions,

- hazards will be avoided,
- workflow will be optimized and accelerated,
- repair costs and machine downtimes will be minimised,
- reliability and durability will increase.

The operating instructions should be read carefully by machine owners and operational officers.

In individual chapters, important safety and hazard instructions are given, by which you are prepared for dangerous situations. It is unfortunately too late to read this manual during operation. In addition, observe the principle:

Careful operation is the best protection against accidents!

The operating instructions are accessible on the work equipment.

2.2. Customer Support

FRANK-PTI feels responsible for all products, which come from our home. Should there be any problem during operation or in the technical area, feel free to contact us at the address listed herein. If you have questions or suggestions for the improvement of our products or require customizations for your particular needs, please con-tact us at:

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Visit our homepage:
<http://www.frank-pti.com> 

2.3. Description

ProbeNet is software for collecting data to the PC from instruments. It is easy to use and displays clearly formatted measurement data. The program is set to receive mode by pressing the start button, and receives the measurement data from all connected devices. Pressing the stop button ends the collection of data.

Clearness:

The data is clearly displayed and is easy to understand. The statistical values for maximum, average, minimum, and standard deviation can all be seen at a single glance. Individual results are listed, and can be excluded from the statistics and reports as anomalies.

Display of curves:

If the results from the measuring devices are returned as curves, these are presented within ProbeNet Mini as diagrams. The display can be enlarged and scrolled to show more detail. This allows study and analysis of interesting areas of the measurement. When new measurements are added, the displayed graphs are automatically adjusted so that all curves are visible.

A hard copy of the report can be printed, to document the results, and contains an area for general information about device and measurement process, a listing of individual results, statistics, and a diagram showing the curves. This area can be removed from the report if desired, in order to save space. A company logo can also be added if required. The results can be exported manually or automatically, both in a spreadsheet (e.g. Microsoft Excel), and a machine-readable format for further processing by other systems. The data can either be saved to the computer's hard disk or saved to a network drive. The program configuration can be password protected, to prevent unauthorised employees making changes to the settings. However all features connected with carrying out the measurements are still available to all users. ProbeNet Mini can be downloaded for evaluation from <http://mini.probenet.biz/>. This is a fully functional version, with support for all the devices of the full version. The only difference is that it can not save data or print reports. Upon purchase of the software you will receive the license key to activate the full version.

2.4. Specifications and technical data

Compatibility:

ProbeNet supports all devices with a data output manufactured and distributed by the PTA Group. If support is not yet available in the program, it can be added if needed. Devices are easily connected, via serial port, USB, or over the network. ProbeNet supports parallel operation of more than one device.

User defined parameters:

To satisfy the demands of various different applications, userdefined parameters can be used to give ProbeNet several profiles. For every test series, the lab worker must select a profile (e.g. customer log, internal log etc.) and data must be entered in the appropriate fields, where it is recorded both as a hard copy report and exported data. To speed data entry, the last values entered are saved, and can be quickly selected if desired. Fields can also be set as optional or required.

System requirements:

Operating System:	Windows XP SP3, Windows 7, 8 8.1 & 10
Software:	Microsoft DotNet 4.0
Resolution:	at least 1280x800 Pixel
Internal memory:	at least 1GB
Hard drive space:	at least 2 GB free space

2.5. Installation


Normally the program is delivered to you saved on a FRANK-PTI USB Stick.

With copy and paste you can just save the whole folder on your desired save path.

Open the ProbeNet.exe file with double-click to start ProbeNet.

2.6. Main menu

After you opened the program the following main menu appears.

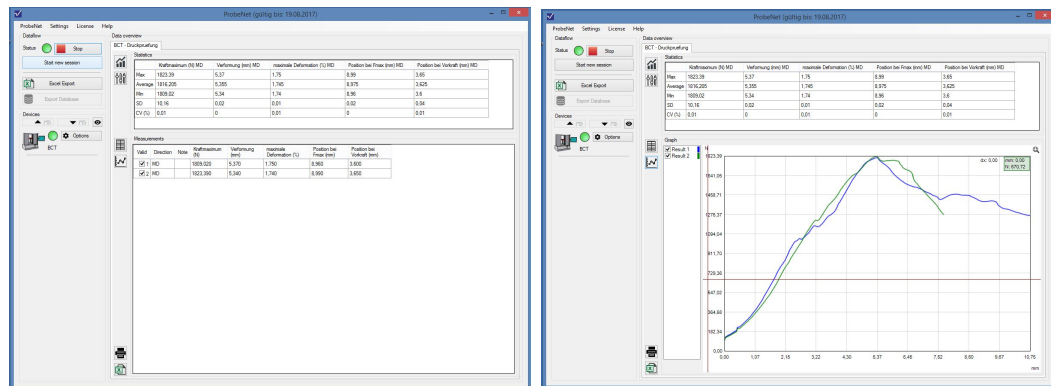
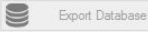
On the left side you can see the rubrics "Data flow" and "Instruments".

Here you can see the status which is indicated by the green or red signal.

With  you can activate the data transmission.

And with  you can finish it.

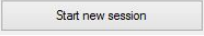
Press  you can generate and start a new measurement series.
The current series will be deleted.

The two buttons below can be used for data output in
excel  or data base .

A window will open which asks you for the destination file path.

For the export to a data base you need to do some pre-settings in the options.

In the rubric instruments the installed and implemented instruments are listed.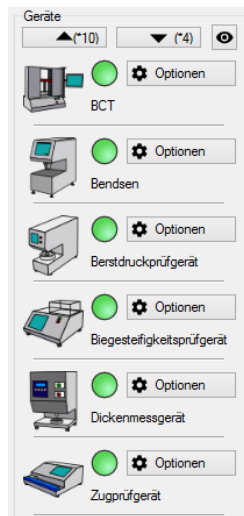

Each instrument is shown with a miniature picture.
The Signal dot in the middle shows the status of each instrument.
If its green this indicates an active data transmission.
Each instrument can be edited with the respective option button.

The arrow buttons can be used to move instruments upwards and downwards to arrange them in your preferred order.
With the eye button you can fade out the single instruments or make them visible again.

On the right, you can see a data overview to a chosen instrument, with individual value sets depending on the machine type.
If the instrument delivers a graph than it is displayed in the lower part.
Also graphs of several measurements can be shown simultaneously in one picture.

You can use the Tabs to switch between the installed instruments and machines.

The upper right window shows the statistics of one series or the used test method and its parameters.


With you can switch to the statistical table view (picture above) and with to the parameters view (picture below).

The lower right window shows an overview of all performed measurement series. (eg. series 1, series 2,...)

With and you can switch between table (picture below) and graphic mode (picture above).

By setting the hooks in the boxes in front of each series single measurement series can be faded out.

By using the mouse you can zoom to single parts the graph. With the magnifier symbol you can zoom out again.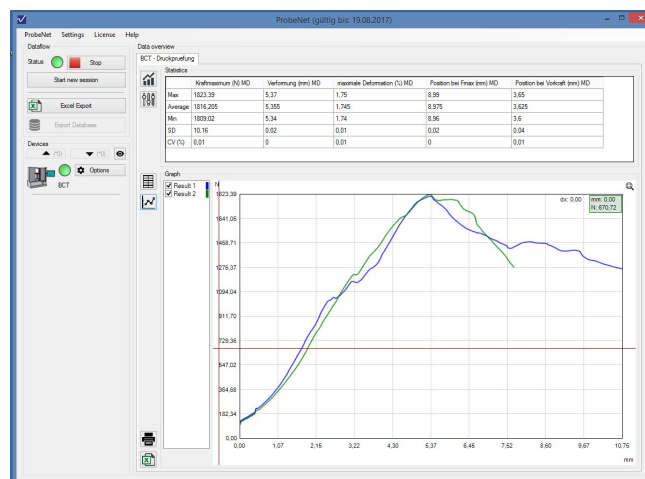


With  you get into the print preview and print settings of the displayed machine where you can print a report.

With  you can export the results of each machine directly into a excel file.


2.7. Menu bar

On the upper end of the main menu you will find the menu bar.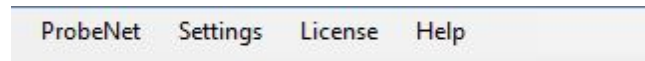

By clicking on the single rubrics a short menu will open.


In the rubric "ProbeNet" you can activate and deactivate the ProbeNet connection with "play" and "hold".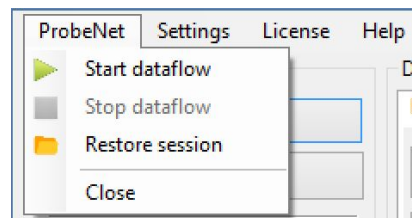


Underneath you can restore a measurement series that was closed before.

With close you can close the sub menu.


In the rubric settings you have the possibility to select the language and to open the settings menu.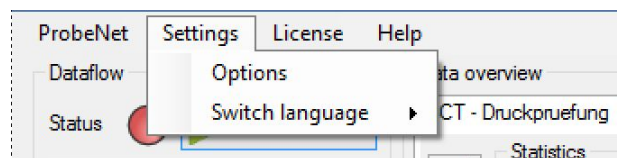

In the rubric "license" you can take an insight to your ProbNet license and also the possibility to manage the devices.

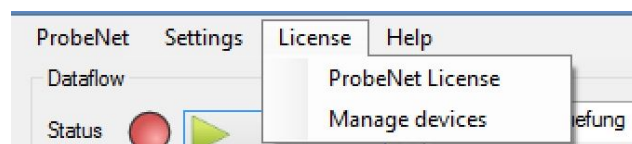
In the rubric "help" you can find a support for troubleshooting, the event log and access to the log folder as well as information about ProbeNet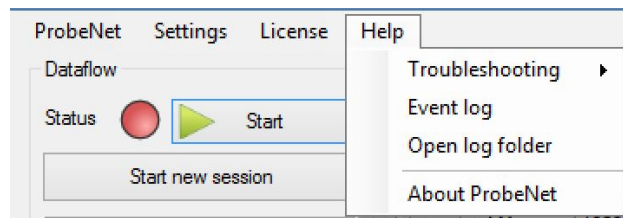


In the trouble shooting sub menu you can check the ProbeNet connection.
When you open "ComPortMonitor" you can edit the baud rate of the serial data transmission and the port.

In the output window the data of the serial transmission will be shown.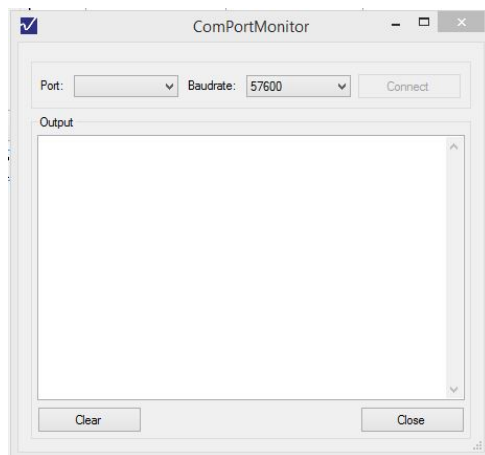

With "clear" you can clear the shown data.
With "close" you will get back to the main menu.

Click on Event log and the following window appears, where you can find the log of the device.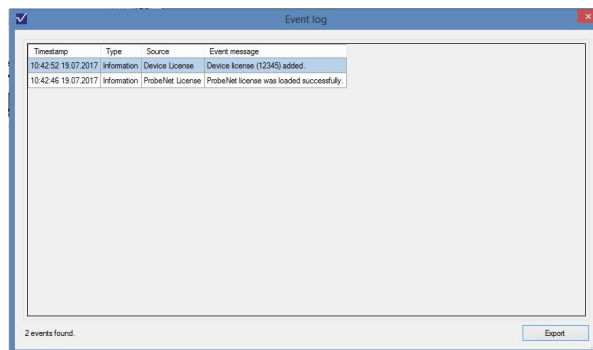


with export you can save the machine log on a desired path.

With "open log folder" you can open the windows folder, in which the logs are saved.


In "About ProbeNet" you can find the installed ProbeNet version and how long your license is valid.
Furthermore you can find the FRANK-PTI Service contact.

2.8. New measurement series

After using the new series button the following window appears.
(optional: only appears if this option is chosen in the settings)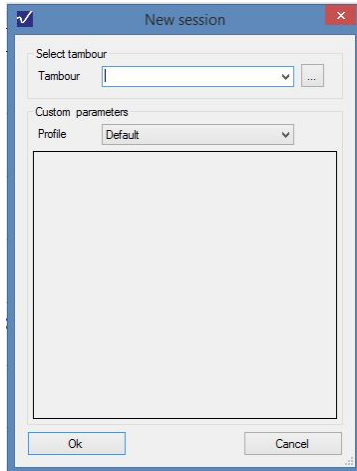

In the upper part you can edit the tambour data.
With the button on the right you get into the tambour settings.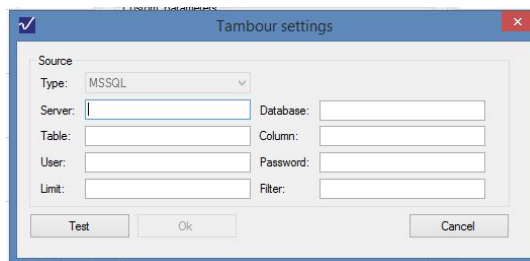

Here you can set and name servers, tables, users limit as well as the data base, rows, passwords and filters.


Now the test button is used to check if the inserted data is correct and can be integrated.

Press OK to confirm.

Press BACK to abort.

2.9. Machine options

Next to each machine in the list you will find a separate option button where you get into the single machines options.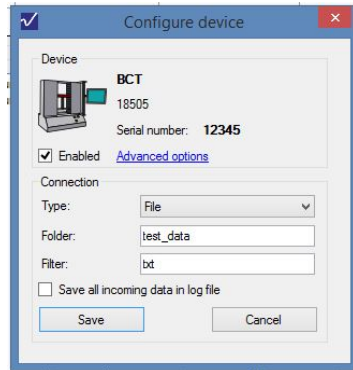


In the upper part you can see the machine data, like serial number and type number.

Here you can edit the connection of each machine.


By setting a hook you can enable the machine.

Press save to fix the currently done changes
Cancel to abort any changes.


Click on the "Advanced Settings" to open the following window: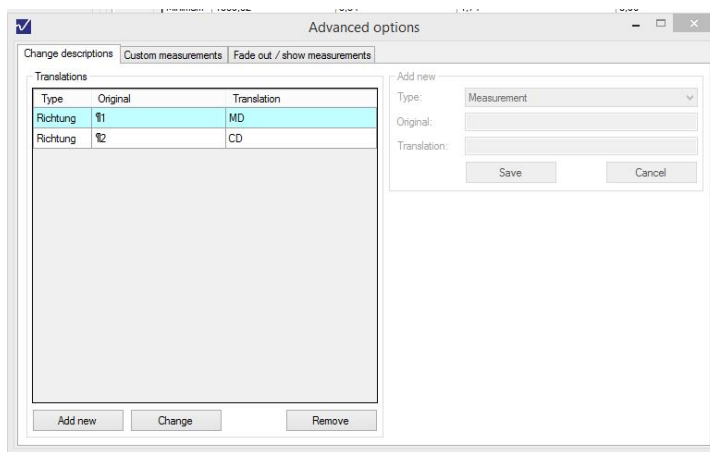

Here you can change descriptions, translate and convert values with a conversion factor.

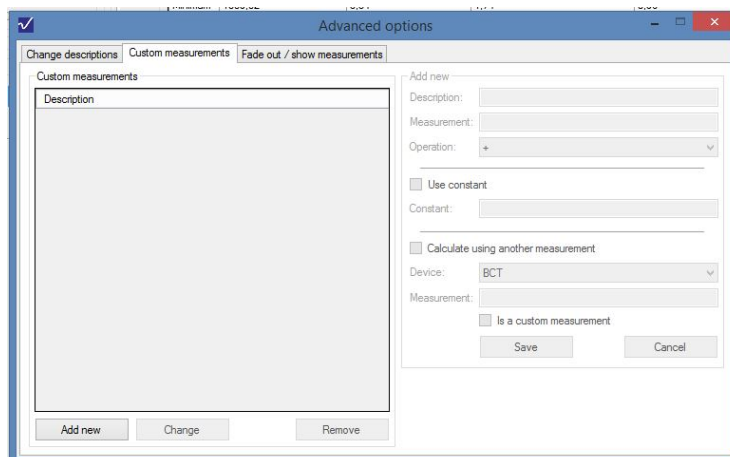
In "Fade out / show measurements" you can determine which measurements and values should be included.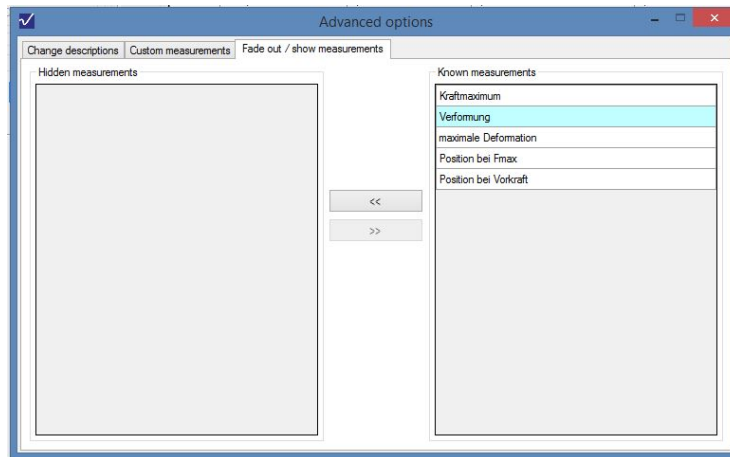


2.10. Options

When you open the settings in the menu bar the following window appears.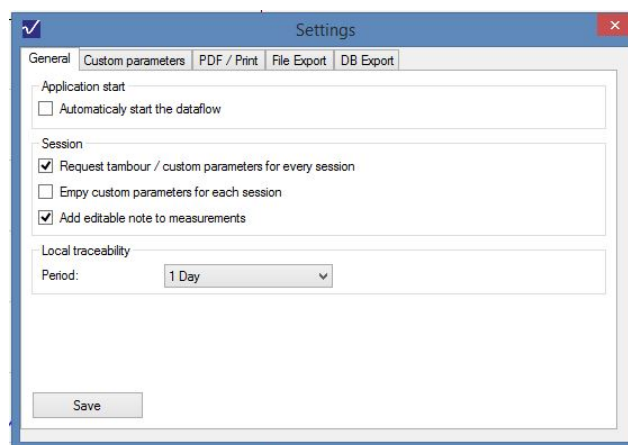


In the tab "general" you can determine if the data transmission starts automatically. You can select if tambour or own parameters should be entered before starting a test series or if the notes should be faded out. Also you can determine over which period of time the data should be saved.

In the tab "custom parameters" you can add new parameters and determine if they are required.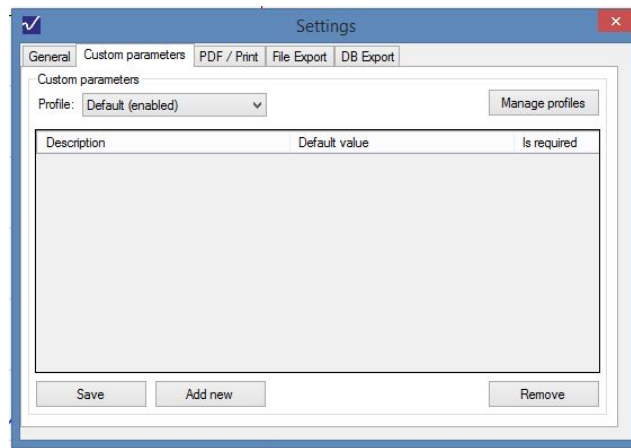


With " add new" you can add a new parameter in the list. And with remove you can remove it.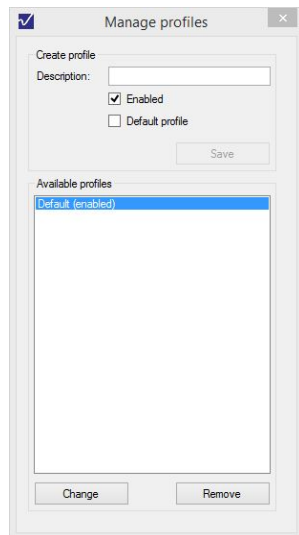

With the button "Manage profiles" you get into the profile menu.


Here you can change and remove profiles.

In the Tab " PDF/print" you can edit the logo that is shown on the printed and PDF- test reports.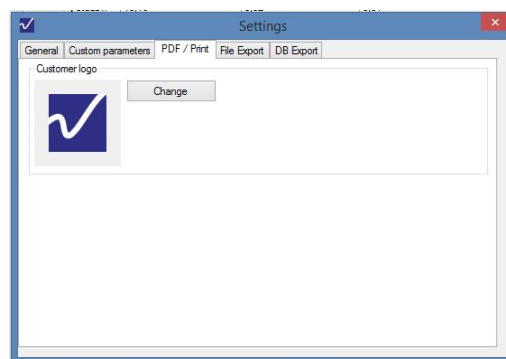


In the "export-tab" you can edit the export path as well as the format of the exported files.

Furthermore you can choose the path of the excel templates.
Press "save" to save the changes.

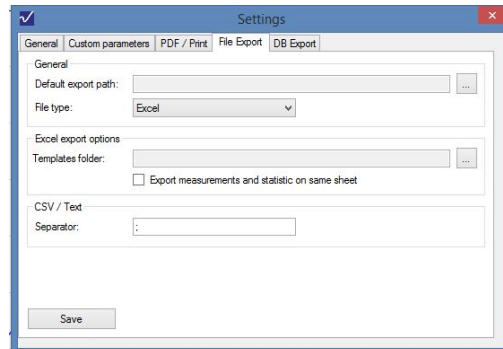
In the tab "DB Export" you can install the connection to a data base set its export options.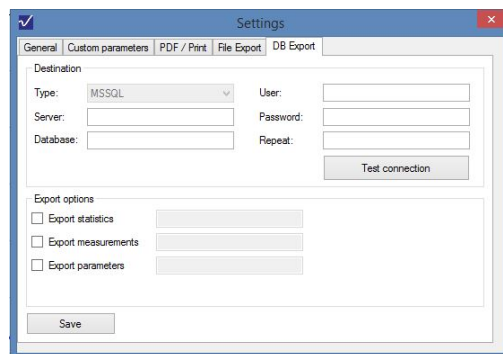

2.11. Print Report

When using the print button the following window opens.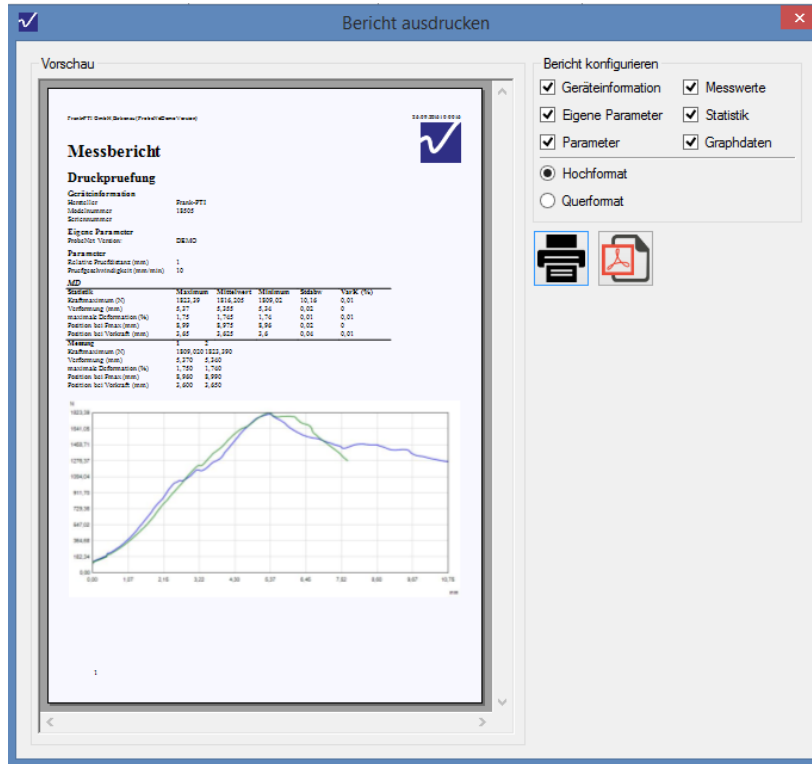

Here you can see a preview of the measurement report.

On the right, you have several options to adapt the report, by setting the hooks for the different rubrics.


Underneath you can choose the format.

With the printer symbol you can print the report directly or export to PDF by using the respective button.

2.12. Create Testsequences for the Modularline


For more information and a detailed description of the ProbeNet software, please refer to the product-specific instructions.

- Start ProbeNet on your PC.
- The main screen appears.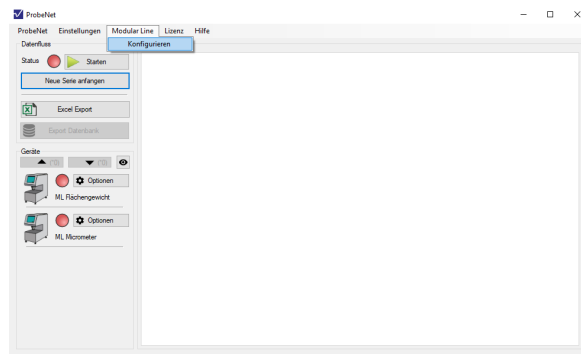


- Click ModularLine in the upper taskbar.
- Click on Configure.

The following page for managing the checking profiles appears: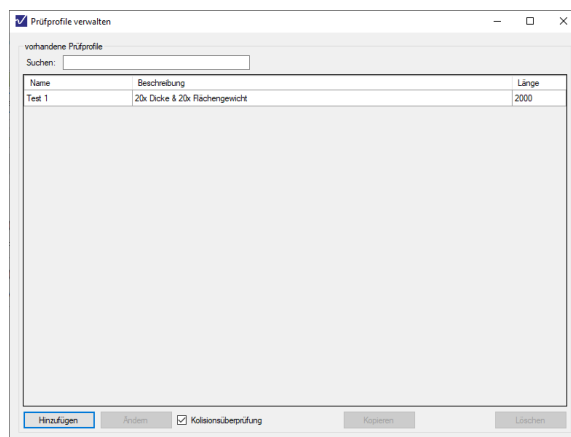


- Click "add" at the bottom left


The following template for creating a new test program appears.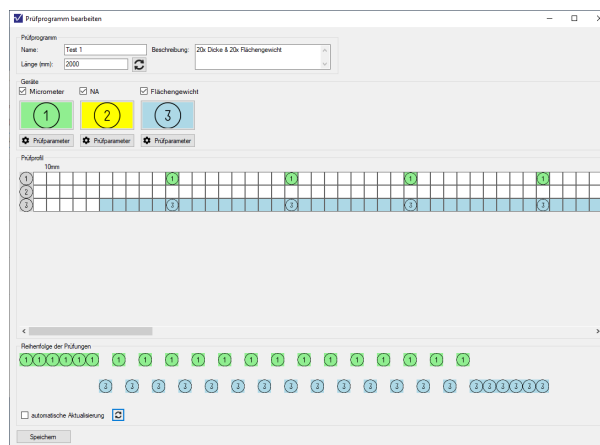


- Set the name, description and length (and click on the "Refresh" button to generate the profile).
- Select the desired test instrument (click on the colored ID)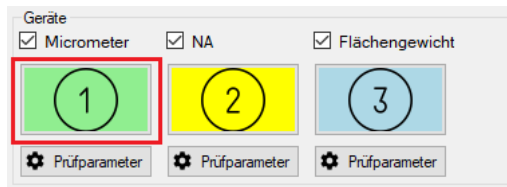


- Click the desired position with the left mouse button.
- Alternatively - with the right mouse button you can set up an automatic placement of the measuring positions: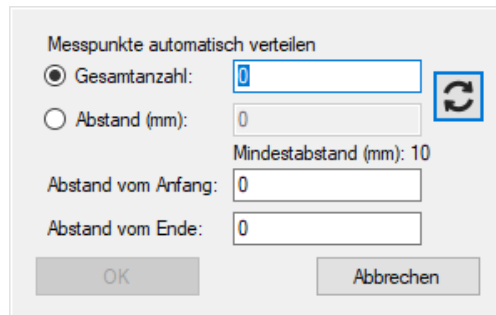

- To remove all measurement positions for the selected device: right-click on the line header: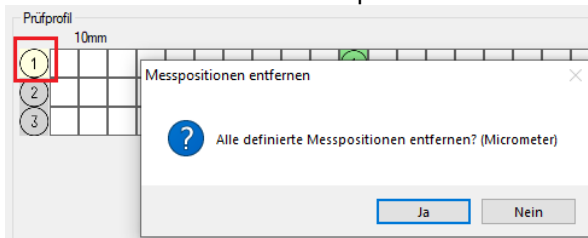

- In the lower area, the chronological order of the tests is displayed:

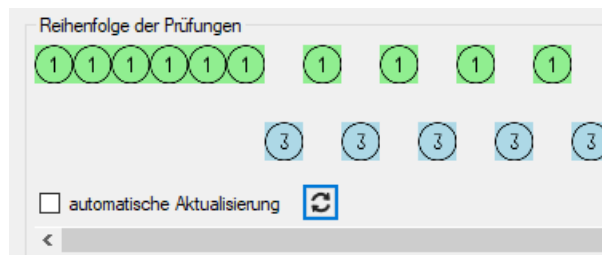- Tests that are carried out at the same time "stand on top of each other".
- I.e. in the area "Test profile" the positions are defined, several test devices could test at the same position. Under "Sequence of tests" you can see the chronological course of the test program.

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All rights reserved. Graphic, electronic or mechanical, including photocopying, recording, taping or the storage and retrieval-systems - without written approval of the editor is prohibited. No part of the document may be reproduced in any form or by any means.

Products that are mentioned in this document are either registered trademarks and/or registered trademarks of the respective owner. The publisher and the author do not assert a claim on these trademarks.

Although this document was prepared with greatest care, errors can be contained in it. In this case we do not take any responsibility for the damages resulting from the use of the document. In no case FRANK-PTI and the author of the document are liable for lost profits or other commercial damages that have supposedly resulted directly or indirectly.

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FRANK-PTI GmbH